

KEOFITT Plastic sleeve for Aseptic Sampling System

ART. NO. 260125

GENERAL



When handling glass bottles, there is always a risk that it may break. The new plastic sleeve offers convenience and also protects the bottle during handling and transport.

The plastic sleeve is a convenient and safe stand/holder for the KEOFITT Aseptic System with 500 ml bottles.

The top part of the plastic sleeve (black) is autoclaved together with the Aseptic System and bottle and after autoclavation the bottom part of the plastic sleeve (transparent) is screwed on to the top part.



It is important that the transparent part of the plastic sleeve is not autoclaved as the material is PVC which will get brittle after autoclavation.

The plastic sleeve is also available for bottles up to 2000 ml, article no 260225.

FEATURES



The top part is attached to the bottle and Aseptic System while the bottom part is screwed on to the top part after autoclavation of the Aseptic System.

CERTIFICATION

• N/A

TECHNICAL DATA


Material (without process contact)

• Top part (black)

POM C

• Bottom part (transparent) PVC



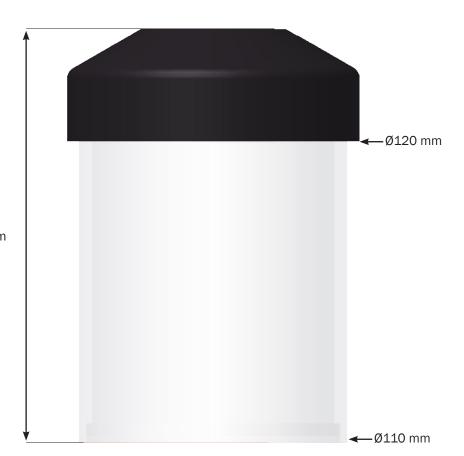

SOME PEOPLE WOULDN'T GAMBLE WITH THEIR SAMPLE™

Data Sheet from keofitt.dk Printed 15 Mar 2010